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(URL <https://www.himac-science.jp>)

Keep this rotor instruction manual and the centrifuge manual in the file.

Swing Rotors for Ultracentrifuge

P55ST2

Before using these rotors, please carefully read this instruction manual, "ROTORS, TUBES, BOTTLES AND CAPS", and the centrifuge instruction manual for its efficient operation and for your safety.

Keep this instruction manual for your reference and refer to it as required.

Contents

1. Specifications	1
2. Applicable tubes/bottles	2
3. How to use	3
4. Maintenance	6
5. Using cesium chloride	8
6. Packing list	9
7. Warranty of the rotor	10
8. Decontamination	12
9. Rotor retirement	12

Part No. { RLM rotor
9122770M
Rotor with optical adapter
91227700
Rotor with optical disk
9122771C

2020.07
99999717

Original instruction manual

Copyright © 2020 Eppendorf Himac Technologies Co., Ltd.

All rights reserved. No part of this document may be reproduced or transmitted in any form or any means without permission from Eppendorf Himac Technologies Co., Ltd.

The names of actual companies and products mentioned herein may be the trademarks of their respective owners.

Safety Reminder

- Centrifuge rotors rotating at high speed have considerable potential for damage to personal properties if used improperly.
For safe and proper use of this rotor, carefully read the centrifuge instruction manual and this rotor instruction manual before use and observe the instructions.

 **WARNING :** and  **CAUTION :** notes are used to call your attention in this manual to prevent personal injury or damage to the rotor and the centrifuge.
These notes are defined as follows.

-  **WARNING:** indicates a potentially hazardous situation which, if not avoided, could result in personal severe injury or possible death.
-  **CAUTION:** indicates a hazardous situation which, if not avoided, could result in personal injury or severe damage to the instrument.

WARNING

- Never use any material capable of producing flammable or explosive vapors. Your centrifuge is not explosion proof centrifuge and your rotor is not explosion proof rotor.
- Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.
- Never exceed the maximum speed of the rotor(mentioned on the rotor).
Always reduce rotor speed as instructed in this manual when rotor speed is limited due to sample density or kinds of tubes.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor(including buckets). Using such a sample could corrode the rotor(including buckets).
- Do not remove the RLM adapter or optical adapter/disk from rotor, or replace it with the adapter/disk for another rotor. The adapter/disk is critical component that detects the over-speed of rotor. If an adapter/disk that is not compatible with the rotor is attached, the rotor could break, resulting in damage to the centrifuge.
- Do not allow the rotor temperature to rise 100°C.: This would cause the material brittle.
- Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.
- When using swing rotor, be sure to set all buckets of the same type whether or not samples are put in buckets: Failure to do so could not only cause the rotor to vibrate, but could result in the rotor being deformed and the buckets becoming detached, which is very dangerous.
- If the centrifuge, rotor, or an accessory is contaminated by samples that toxic or radioactive, or blood samples that pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the centrifuge, rotor, or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the centrifuge, rotor, or the accessory properly before requesting repairs from an authorized sales or service representative.
- It is your responsibility to sterilize and/or decontaminate the centrifuge, rotor, or parts properly before returning them to an authorized sales or service representative.

CAUTION

- Check the chemical resistance chart attached to the rotor, and do not use any sample inapplicable to the tubes or the tube adapter. Using such a sample could deteriorate them.
- The allowable speed may be lower depending on the density of sample, use of salt-such as cesium chloride-or the combination of tubes and tube adapters. Refer to this instruction manual.
- Always balance the tubes and samples within the allowable imbalance of rotor, including the tube adapters. Do not exceed the allowable imbalance. Otherwise the rotor and the centrifuge may be damaged.
- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.
- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Use the rotor tubes and bottles within their actual capacities.
- Do not use tubes that have exceeded their life expectancy. Failure to do so could result in damage of tubes and the rotor and the centrifuge. The life expectancy of tubes depends on factors such as the characteristics of samples, speed of the rotor used, and temperature. Always check for deterioration and damage (cracks, deformation, and so on) on tubes before using them. Do not use the tubes if you find such a problem.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
Contact an authorized sales or service representative.

1. Specifications

1.1 Construction

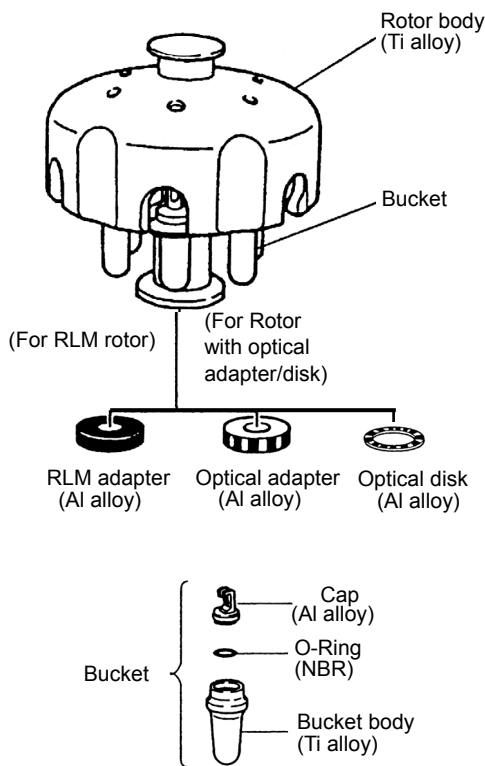
Fig.1-1. Construction of P55ST2 rotor

1.2 Specifications

- Max. speed: 55,000 rpm
- Max. RCF: $366,000 \times g$
- Allowable density at maximum speed of the rotor: 1.2g/mL
- K Factor at maximum speed of the rotor: 50
- Nominal rotor capacity: 5 mL \times 6 tubes = 30 mL
- Material of rotor: Ti alloy
- Tube size: $\phi 1.3 \times 5.2$ cm
- Weight: 5.5 kg
- Max. diameter: 16.5 (21.8) cm
- Max. height: 18.2 cm

(The numerical value in parentheses is maximum. diameter measured with the bucket swinging.)

⚠ CAUTION:

- Do not use a RLM rotor with the centrifuge other than a series of a CP-NX/WX/MX/ α ultracentrifuge.
- Do not use rotor with optical adapter/disk with model 55P and 40P (old type centrifuge).

1.3 Cross-sectional view of rotor

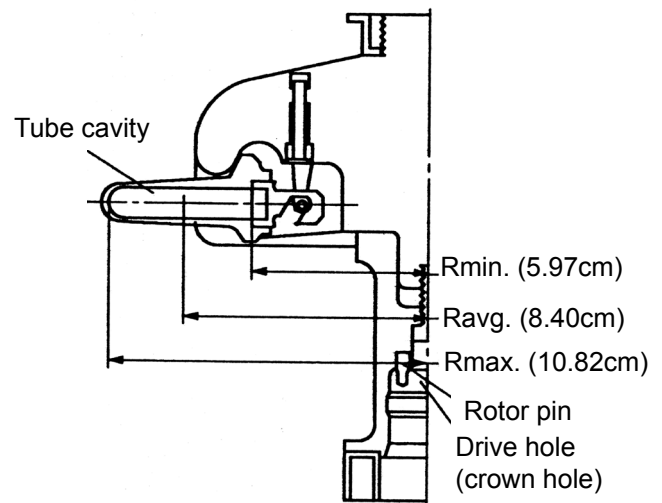
Fig.1-2. Cross-sectional view of P55ST2 rotor

⚠ CAUTION:

Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

1.4 Characteristic of rotor

Table 1-1. Characteristic of P55ST2

Rotor Speed (rpm)	RCF ($\times g$)			K factor
	Rmax.	Ravg.	Rmin.	
10,000	12,100	9,390	6,670	1,506
20,000	48,400	37,600	26,700	377
30,000	109,000	84,500	60,100	167
40,000	194,000	150,000	107,000	94
50,000	302,000	235,000	167,000	60
55,000	366,000	284,000	202,000	50

※For detail about K factor, refer to “ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)” attached to the centrifuge.

Calculation formula

$$RCF = 1.118 \times 10^{-5} \times R \times N^2 (\times g)$$

R: Rotating radius (cm)

N: Speed (rpm)

2. Applicable tubes/bottles

2.1 Applicable tubes/bottles

Table 2-1. Applicable tubes/bottles

Nominal capacity (mL)	Tubes/Bottles				Tube adapter		Cap		Max. speed (rpm)	Max. RCF (xg)
	Part No.	Name	Size (φ X L cm)	Actual capacity (mL/tube)	Part No.	Name	Part No.	Name		
5	332245A	5PA tube	1.3 X 5.2	4.9	---	---	---	---	55,000	366,000
	S303273A	5PET tube	1.3 X 5.2	5.3	---	---	---	---	55,000	366,000
3	S301599A	3PC tube	1.3 X 5.1	3.4	---	---	---	---	55,000	366,000
0.9	S304296A	0.9PC thick-walled tube	0.8 X 5.0	0.7	336697A	2S5 adapter	---	---	40,000	189,000

* The material of tube adapter: Polyacetal


- ⚠ CAUTION:**
- When using 2S5 adapters with 0.9PC thick-walled tubes, run the rotor at 10°C or less, otherwise the tubes deform during operation. As a result, it can be impossible to remove the tubes from 2S5 adapters.
 - When using 0.9PC thick-walled tubes with 2S5 adapters, 0.9PC thick-walled tubes are disposable.
 - Chemical resistance and the strength of tubes varies with speed, temperature, and so on. Before using sample, fill tubes with water, buffer solution, etc. instead of sample and run them at the intended speed, temperature, and so on to ensure that there is no abnormality.

Refer to “ROTORS, TUBES, BOTTLES AND CAPS (Part No. S999204)” for the cleaning and sterilization of the tubes/bottles.

3. How to use

How to use tubes/bottles varies in the kinds of tubes/bottles. The brief procedure of tubes/bottles handling is described as follows.

For the detail about handling tubes/bottles, refer to “ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)” attached to the centrifuge.

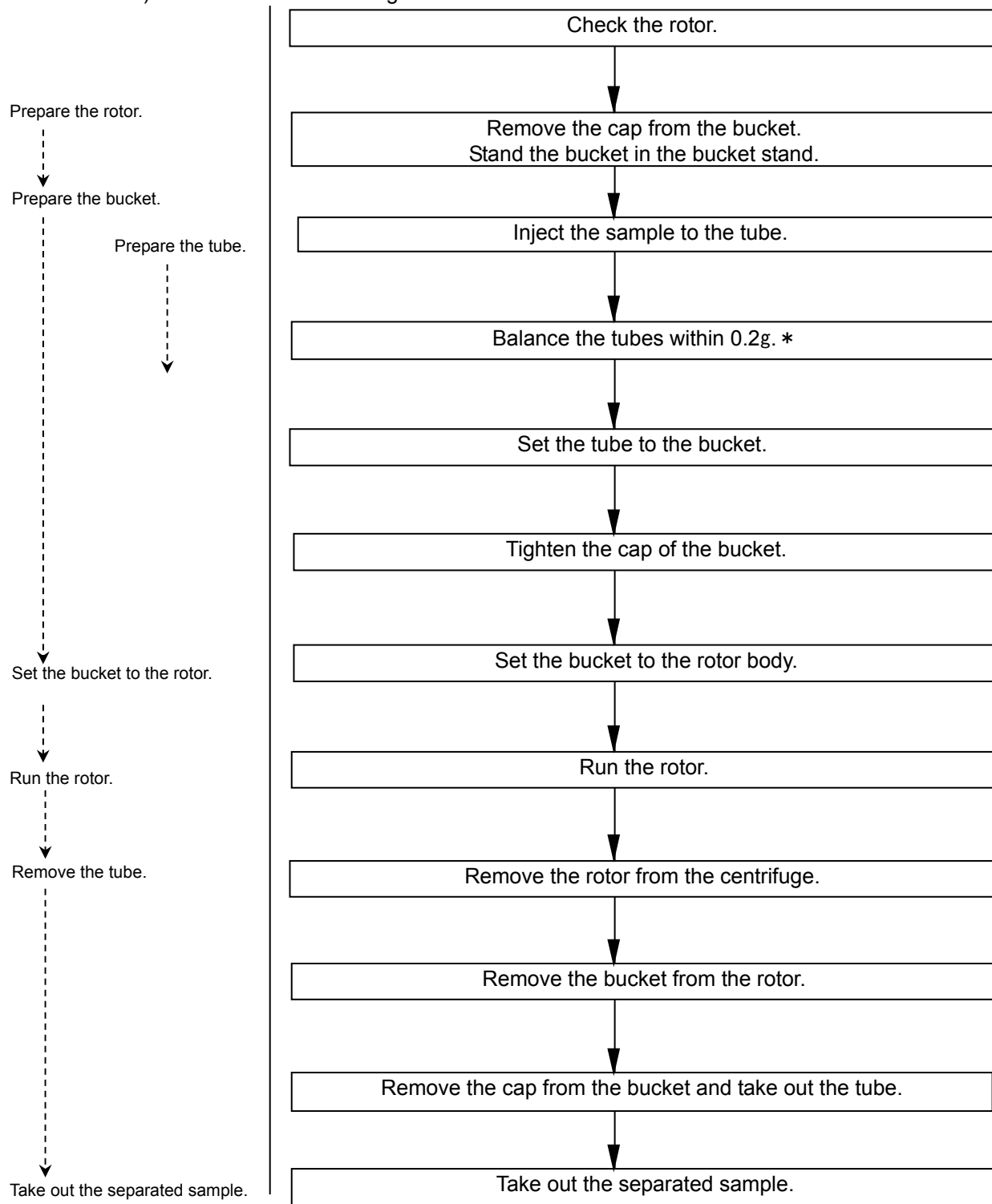(*) When using CP-NX series, CP-WX series, CP-MX series, or CP- α/β series, approximately even quantities of sample in the tubes are sufficient for balancing, and extremely different sample quantities must be avoided. (The levels of samples in the tubes should be approximately equal. Their difference should be within 5mm.)

For some kinds of tubes, if a small quantity of a sample is injected to them, the rotor can not be used at the allowable maximum speed. Tube deformation may occur due to a small quantity of a sample during operation.

=For handling=

● Allowable speed of rotors

1.To centrifuge a sample of an average density more than 1.2g/mL.

Determine the allowable speed from the following equation.

$$\text{Allowable speed(rpm)} = 55,000(\text{rpm}) \times \sqrt{\frac{1.2 (\text{g} / \text{mL})}{\text{Mean density of sample(g/mL)}}}$$

2. The allowable speed may be lower depending on the density of a sample when using of salt-such as cesium chloride. Refer to “5. Using cesium chloride” for the detail.

3. The allowable speed may be lower depending on the combination of tubes and tube adapter, etc. Refer to “2. Applicable tubes/bottles”

4.Reduce the rotor speed when the rotor has reached the primary life. Refer to “4. Maintenance●Life of rotor”.

●Tube/Bottle

Before operation, check whether the combination of tubes and tube adapter is properly or not.

Fill the tube up to within 3mm from the top of the tube with sample.

Refer to “2. Applicable tubes/bottles” in this manual and “ROTORS, TUBES, BOTTLES AND CAPS(Part No. S999204) ”.

●Handling the rotor

WARNING :

●Never exceed the maximum speed of the rotor(mentioned on the rotor). Otherwise, rotor may be damaged and result in damage to the centrifuge.

●Check the chemical resistance chart attached the rotor (including buckets), and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor(including buckets).

CAUTION :

●Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.

●Balance the tubes/bottles and the samples including the caps and the tube adapters within the allowable imbalance of the rotor.

Do not exceed the allowable imbalance. Otherwise, the centrifuge or the rotor may be damaged due to excessive imbalance operation.

●If foreign matter adheres to the rotor(including buckets), wipe it off promptly. Otherwise, the rotor(including buckets) may be corroded.


Set the balanced tubes to the rotor symmetrically.

<Tightening the cap to the bucket>

CAUTION :


Imperfect tightening of the cap may cause imbalance operation and it may happen the serious accident.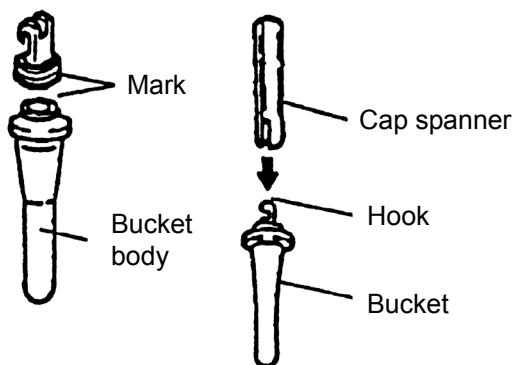

Tighten the numbered cap to the same numbered bucket with the supplied cap spanner to match each mark of the bucket and the cap.

Tighten the cap until the bottom of the cap bumps to the bucket and the cap must be adjusted to the match mark of the bucket.

<Setting the bucket to the rotor>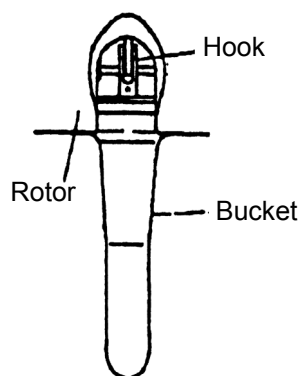

Hang the hook of the bucket on the pin of the rotor.


Match the numbered buckets with number of rotor and set the buckets to the rotor.

WARNING :

Be sure to set all buckets of the same type whether or not samples are put in buckets: Failure to do so could not only cause the rotor to vibrate, but could result in the rotor being deformed and the buckets becoming detached, which is very dangerous. Never use buckets made by other companies, or any other type of bucket that is not exclusively made for the rotor even if it is made by us.

< Examples of incorrect bucket setting >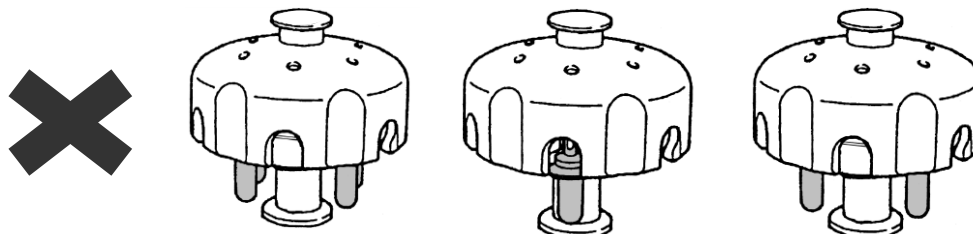

Tighten the cap to the bucket securely. Set the bucket to the rotor securely.

WARNING :

Before setting a rotor in ultracentrifuge, make sure that buckets have been securely installed in the rotor: Incorrect installation of buckets in rotor could damage the rotor or detach the buckets, which is very dangerous.

CAUTION :

Inspect and maintain the rotor after use. If abnormality is observed, do not use it. Contact an authorized sales or service representative

● Sample limitation

WARNING :

- Never use any material capable of producing flammable or explosive vapors. Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage.
- When using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.

● The rotor life management of the rotor with optical adapter/disk

CAUTION :

Each time the rotor with optical adapter/disk is used, be sure to enter the results in the “rotor log book”. Failure to do so will disable management of the life of rotor.
No entry is needed for RLM rotor.

● RLM adapter and Optical adapter/disk

WARNING :

Do not remove the RLM adapter or optical adapter/disk from rotor, or replace it with the adapter/disk for another rotor. The adapter/disk is a critical component that detects the over-speed of rotor : If an adapter/disk that is compatible with the rotor is attached, the rotor could break, resulting in damage to the ultracentrifuge.

CAUTION :


- Do not bring the RLM adapter near a magnet, or scratch it: Doing so will erase the memory stored in adapter, and make the rotor unusable. To prevent the adapter from being scratched, store the rotor with RLM adapter, using the stand provided with the rotor (rotor stand for protecting adapter). If the adapter is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- Take great care not to scratch the optical adapter/disk: Any scratch on adapter/disk will make the rotor unusable. To prevent the adapter/disk from being scratched, store the rotor with optical adapter/disk, using the stand provided with the rotor (rotor stand for protecting adapter/disk). If the adapter/disk is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- If sample etc. adheres to the RLM adapter or optical adapter of rotor, immediately wipe it off with a soft cloth etc. If the RLM adapter or optical adapter corroded, the rotor will be unusable.

4. Maintenance

●Maintenance of rotor

⚠ WARNING :
Do not allow the rotor temperature to rise over 100°C.
Otherwise, the rotor can be brittle.

⚠ CAUTION :
Use a neutral detergent having a pH between 5 and 9.
Otherwise, the rotor can be corroded.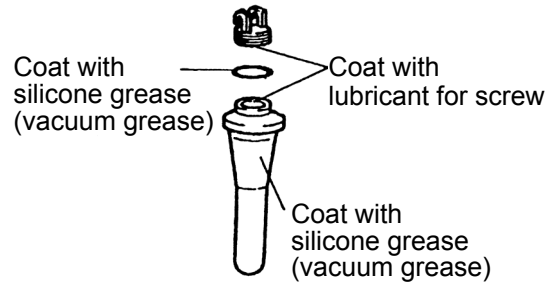

Wash the bucket with tap water or a dilute solution of neutral detergent and rinse it out with distilled water. Turn the bucket upside down with the cap removed to dry it well. Check the bucket is completely dried and then put the light coat of silicone grease (vacuum grease) on the surface of the bucket, the light coat of lubricant for screw on the thread portion, the light coat of silicone grease (vacuum grease) on the O-ring and the gasket. Store them in a dry place. If foreign matter adheres to the bucket, soak the bucket in warm water in one or two hour and remove foreign matter with a soft brush etc. Use a neutral detergent having a pH between 5 and 9. If you can not remove it, contact an authorized sales or service representative. Wipe the stain off the rotor body and then put the light coat of silicone grease (vacuum grease) on the surface of the rotor body. Store it in a dry place. Put the light coat of silicone grease (vacuum grease) inside the drive hole(crown hole) and then wipe it out with a soft cloth.

●Sterilizing rotor

⚠ WARNING :
Never sterilize the rotor(including buckets) by autoclaving or boiling. Otherwise the strength of rotor(including buckets) may be significantly decreased.

⚠ CAUTION :
Do not sterilize the tube adapters made of plastic by autoclaving or boiling. Otherwise the tube adapters may be deformed. Such the tube adapters can not be used.

Sterilize the rotor according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin).

Sterilize the tube adapters made of plastic according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin).

* Do not dip the rotor and the tube adapters in the formalin (3%) solution more than 2 hours.

●Inspecting rotor

Check the rotor after every run as follows.

- The thread portion is worn. Contact an authorized sales or service representative.
- The tube cavity is corroded. Contact an authorized sales or service representative.
- The O-ring, the gasket, etc. are deteriorated. Replace it with the new one.
- The adapter is stained. Wipe it out with a soft cloth.

Check the rotor every 100 hours as follows.

- The rotor pin is bent. Contact an authorized sales or service representative.

Check the rotor once a month as follows..

- The inside of drive hole(crown hole) is stained. Wipe it out with a soft cloth dampened with water.

⚠ CAUTION :
Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

●Life of rotor

While using rotor repeatedly, it's strength decreases gradually due to fatigue and creep of material, by which the rotor life is determined.

Sum up the number of operations and number of operating hours. And if either value reaches the figures (the primary life) shown in Table 4, decrease the maximum speed by 10%.When the rotor has reached the primary life, maintenance of the rotor is required by us (extra-cost option).If the operations or operating hours of rotor reach the figures in Table 4 after decreasing the maximum speed by 10%(Secondary life),don't use the rotor any more. If you use the RLM rotor at a speed lower than its maximum permissible speed, the life of the RLM rotor will be prolonged automatically by calculating from the strength of the rotor material.

Read "7.Warranty of the rotor" carefully.


Table 4-1. Rotor life

1,000 operations	2,500 hr
------------------	----------

● Replacing the optical disk

(1) Preparation

- Prepare a new optical disk and knife.
Make sure that the number of stripes of the new optical disk is right as follows.
- Wash and then dry the rotor well.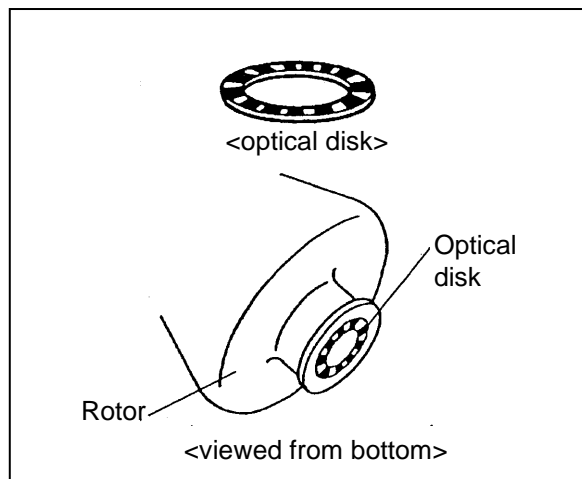

Rotor's maximum speed and optical disk's number of stripes

Rotor's maximum speed (rpm)	Optical disk's number of black bands	Rotor's maximum speed (rpm)	Optical disk's number of black bands	Rotor's maximum speed (rpm)	Optical disk's number of black bands
100,000	9	50,000	18	30,000	29
90,000	10	48,000	19	28,000	32
80,000~83,000	11	45,000	20	27,000	33
70,000	13	42,000	21	25,000	35
65,000	14	40,000	22	23,000	39
60,000	15	38,000	24	21,000	40
55,000~56,000	16	35,000	25	19,000	45
54,000	17	32,000	28	17,000	50

(2) Hold the rotor upside down.

(3) Pry the edge of the optical disk with the knife and remove the disk. Be very careful not to damage the rotor.

(4) Clean the disk hole on the rotor by using alcohol.

(5) Remove the backing paper from the new optical disk.

Place the optical disk so that it snugly fits into the groove of the disk hole on the rotor. Make sure the disk does not move.

5.Using cesium chloride

The cesium chloride(CsCl) solution is frequently used as a density gradient medium, but if the CsCl solution with a high density is used, it may saturate during the rotation of the rotor depending on the rotation condition which may cause CsCl crystals to be educed. Educated crystals have a high density(approx.4g/mL) and apply an excessive load to the rotor which is very dangerous.

The crystallization will also greatly affect the density gradient of the CsCl solution and separation status of the sample. Therefore, always use a solution which will not crystallize the allowable speed.

Fig .5-1 shows the relationship between the maximum allowable density and the speed at which cesium chloride does not crystallize in the condition that standard tubes is filled with the cesium chloride solution. In this diagram, crystallization will occur above the curves.

Be sure to select the density of the cesium chloride solution and the rotor speed at which cesium chloride does not crystallize. The curve in this diagram vary in temperature etc. Before actual operation, centrifuge the cesium chloride solution with the same density to ensure that crystallization will not occur.

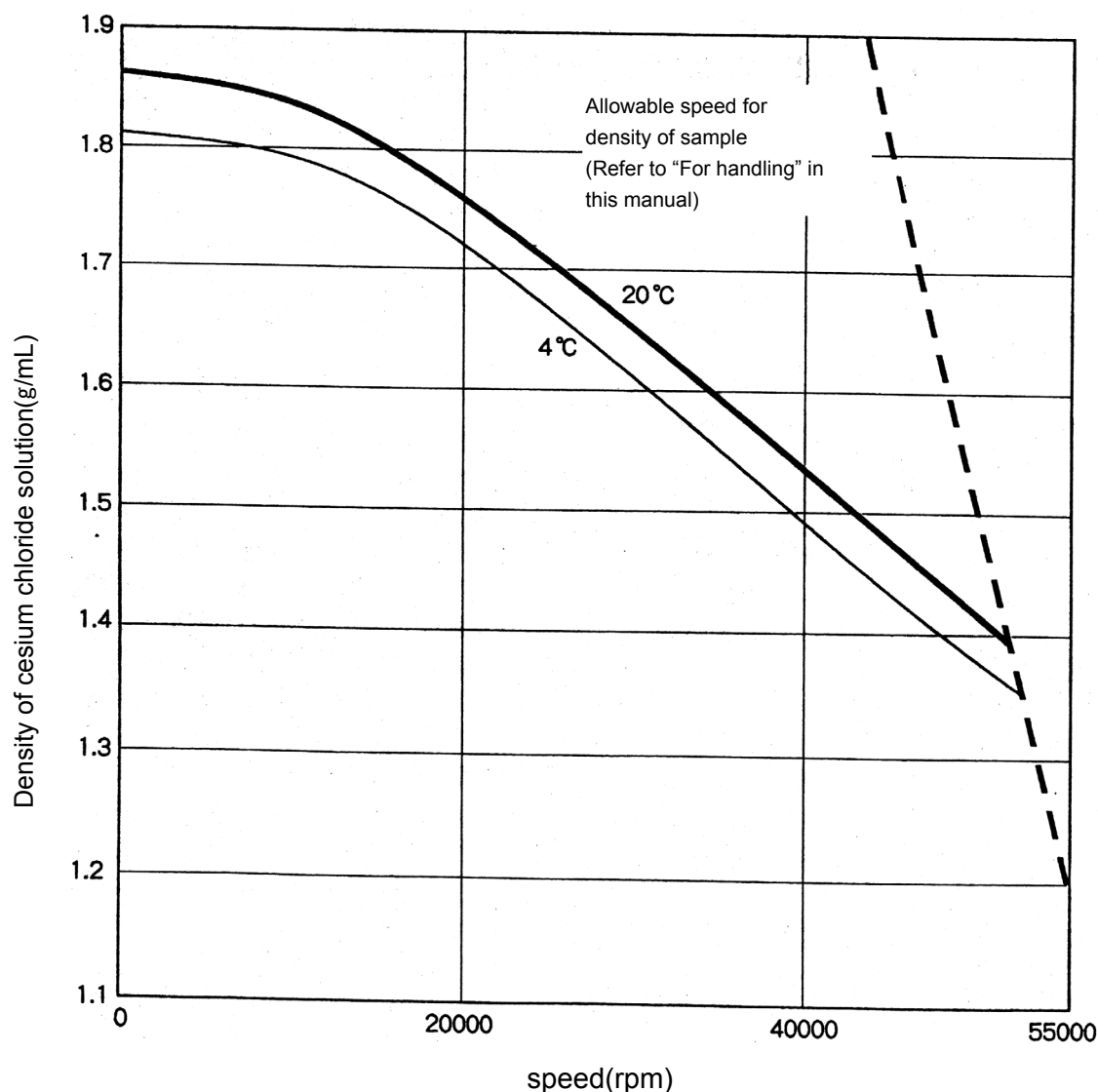
Fig.5-1. Relationship between density of cesium chloride and speed

6. Packing list


Check the parts and accessories of rotor when receiving the rotor.
Contact an authorized sales or service representative when the need arises.

No.	Name	Qty	Remark	Parts No. (Qty)	No.	Name	Qty	Parts No. (Qty)
①	Rotor body	1			⑪	Rotor stand	1	216715 (1)
②					⑫	Lubricant for screw	1	84810601 (1)
③	Optical disk *2	2	For spare	3464445	⑬	Silicone grease (vacuum grease)	1	483719 (1)
④					⑭	Tool box(C)	1	
⑤	Bucket	6			⑮	Cap spanner	1	S407836 (1)
⑥	Bucket body	6*1			⑯	Tweezers	1	452701 (1)
⑦	Cap	6*1			⑰	Bucket stand	1	208204B (1)
⑧	O-Ring	6*1		S401822A (12)	⑱	Rotor log book *3	1	999494 (1)
⑨	O-Ring	18	For spare		⑲	Instruction manual	1	999997 (1)
⑩	5PA tube	2	50pcs./set	332245A(50)	⑳			

*1: the quantities of the parts that are mounted to the rotor

*2: The "optical disk" is only provided with the rotor destined for Europe and U.S.A.

*3: The "rotor log book" is not included in RLM rotor.

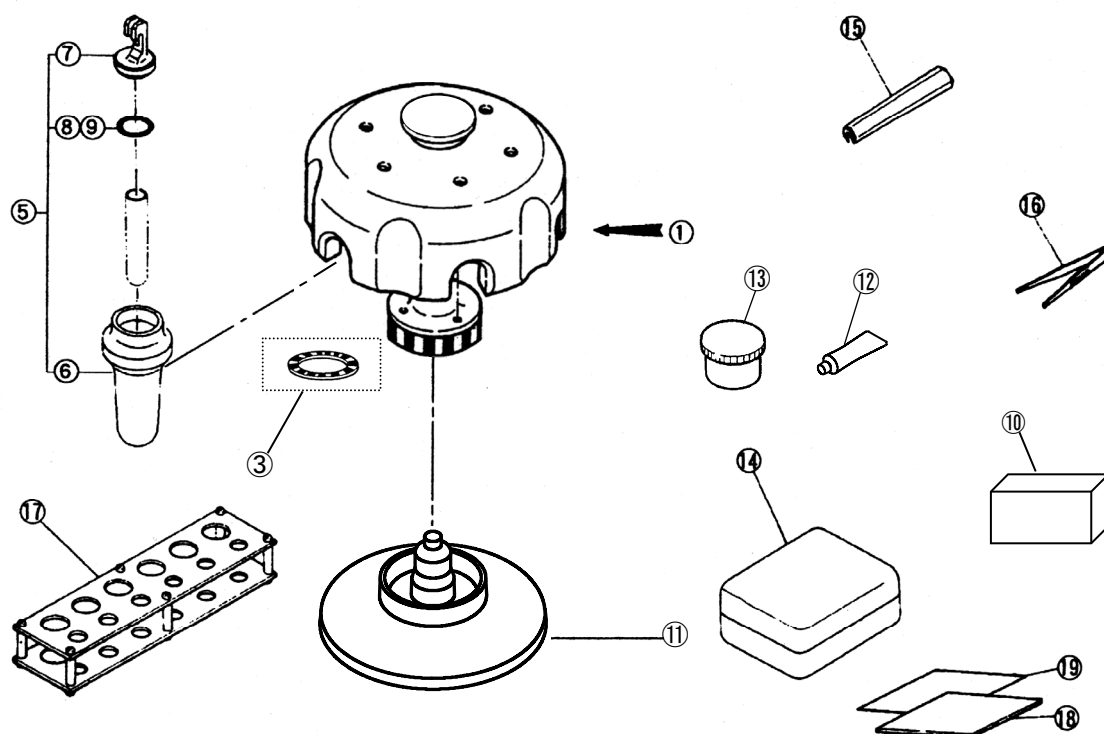7. Warranty of the Rotor

●Warranty for Ultracentrifuge Rotors

Eppendorf Himaс Technologies warrants its ultracentrifuge rotors under the terms and conditions set forth below.

1. Scope of Application

This warranty applies to:

- (1) failure or damage to the rotor due to rotor defects, in materials or workmanship, that are attributable to Eppendorf Himaс Technologies, and
- (2) failure or damage to an Eppendorf Himaс Technologies ultracentrifuge due to defects in a rotor covered by the warranty.

Note: This warranty covers the rotor or ultracentrifuge only, and Eppendorf Himaс Technologies shall not be liable for damage to samples or other items that may arise from failure of or damage to the rotor or ultracentrifuge.

2. Warranty Period

This warranty is valid for five years from the date of delivery of the rotor by Eppendorf Himaс Technologies or an Eppendorf Himaс Technologies dealer. However, except as provided under (1) ②below, even within such period the warranty does not cover cases where the user has exceeded the applicable limit on the number of runs or the cumulative operating time for the type of rotor, as set forth in (1) and (2) below.

(1) Aluminum alloy angle rotor, swing rotor, zonal rotor, and continuous flow rotor

- ① The limits on the maximum number of runs and the maximum cumulative operating time for these rotor types are stated in the table below. The warranty covers the time up until either of these limits is reached, provided that it is reached during the warranty period and that the maximum and allowable speeds are not exceeded.
- ② In order for the warranty to cover use after either of the limits in the table below is reached, inspection by Eppendorf Himaс Technologies will be necessary. The user must pay for such inspection.

If Eppendorf Himaс Technologies' inspection finds that the rotor is free of corrosion and other abnormality, Eppendorf Himaс Technologies will then further warrant the rotor up until either of the limits in the table below is reached once more, but with the rotor's maximum speed reduced by 10%.

Even in such case however, the warranty period shall remain limited to five years from the date of delivery and shall not be extended.

Maximum number of runs	Maximum cumulative operating time
1,000	2,500 hours

(2) Titanium alloy angle rotor, titanium alloy vertical rotor, and neo-angle rotor

The limits on the maximum number of runs and the maximum cumulative operating time for these rotor types are stated in the table below. The warranty covers the time up until either of these limits is reached, provided that it is reached during the warranty period and that the maximum and allowable speeds are not exceeded.

(With these types, the warranty cannot be extended beyond the stated limits, even if the rotor undergoes inspection by Eppendorf Himaс Technologies.)

Maximum number of runs	Maximum cumulative operating time
5,000	10,000 hours

3. Warranty Terms

- (1) Should a rotor covered by this warranty fail or be damaged, Eppendorf Himaс Technologies will replace it with a new rotor in return for payment of a sum determined according to the formula below, based on the ratio of the actual number of runs or cumulative operating time to the maximum number of runs or cumulative operating time.

$$[\text{Replacement price}] = [\text{Current price of rotor}] \times \left[\frac{\text{Actual number of runs}}{\text{Max. number of runs}} \text{ or } \frac{\text{Actual cumulative operating time}}{\text{Max. cumulative operating time}}, \text{ whichever is greater} \right]$$

- (2) Should an ultracentrifuge fail or be damaged due to failure of or damage to a rotor covered by this warranty, Eppendorf Himaс Technologies will supply, free of charge, all parts required for repair of the ultracentrifuge.

However, in the case of failure of or damage to an ultracentrifuge's drive unit, Eppendorf Himaс Technologies will replace it with a new drive unit in return for payment of a sum derived by multiplying the current price of the drive unit by the ratio of the actual total number of revolutions or years completed to the warranted total number of revolutions or warranted years that are stated in the operator's manual for the ultracentrifuge, as in the formula below.

$$[\text{Replacement price}] = [\text{Current price of drive unit}] \times \left[\frac{\text{Actual number of revolutions}}{\text{Warranted total number of revolutions}} \text{ or } \frac{\text{Years completed}}{\text{Warranted years}}, \text{ whichever is greater} \right]$$

4. Notes

- (1) This warranty extends only to the original buyer from Eppendorf Himaс Technologies or the original buyer from an Eppendorf Himaс Technologies dealer. Rotors whose ownership passes from the original buyer to a third person due to resale, transfer or the like are not covered by this warranty. Neither does this warranty cover cases where the ownership of a rotor is reacquired subsequently by the original buyer.
- (2) The warranty for the rotor proper (warranty period and terms) shall continue to apply in cases where rotor parts and the like are additionally purchased.
In such cases, the warranty period shall continue to be measured from the date of delivery of the rotor. If such parts are subject to a restricted warranty period, this warranty shall only cover failure or damage involving such parts that occurs both within the rotor warranty period and within the warranty period for the parts concerned.
- (3) This warranty does not cover any of the following, even during the warranty period:
- ① Failure or damage due to abuse or misuse.
 - ② Failure or damage due to operation or maintenance in a manner contrary to the instructions in the operator's manuals for the rotor and ultracentrifuge.
 - ③ Failure or damage due to disassembly or modification without the permission of Eppendorf Himaс Technologies.
 - ④ Failure or damage due to use of a bucket, adapter, tube, bottle, or the like, other than those approved by Eppendorf Himaс Technologies for use with the product.
 - ⑤ Failure or damage due to disaster such as fire or earthquake.
 - ⑥ Failure or damage due to use with an ultracentrifuge not of Eppendorf Himaс Technologies manufacture without the permission of Eppendorf Himaс Technologies.
 - ⑦ Failure or damage due to use with an ultracentrifuge that has been modified without the permission of Eppendorf Himaс Technologies.
 - ⑧ Failure or damage due to a failed or damaged part (including a tube, bottle, adapter, etc.) not submitted to Eppendorf Himaс Technologies.
 - ⑨ Failure or damage not recorded in a rotor use log book (in the case of an optical adapter type rotor).
 - ⑩ Replacement of consumable items.
 - ⑪ Parts subject to a restricted warranty period that has expired.

●Other Matters

Rotors that are delivered to Eppendorf Himaс Technologies for inspection or repair must be sterilized and cleaned beforehand.

Be aware that Eppendorf Himaс Technologies will sometimes decline to carry out inspection or repair if a rotor or other related equipment is or could be contaminated as a result of use without protection or under inadequate safety management in a radioisotope facility or a laboratory of the level P2 or higher.

8. Decontamination

WARNING:

- If the centrifuge, rotor or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the rotor or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the rotor or the accessory properly before requesting repairs from an authorized sales or service representative. Note that we cannot repair the centrifuge, rotor or the accessory unless sterilization or decontamination is completed.
- It is your responsibility to sterilize and/or decontaminate the rotor or parts properly before returning them to an authorized sales or service representative. In such cases, copy the attached decontamination sheet and fill out the copied sheet, then attach it to the item to be returned. We may ask you about the treatment for the rotor or parts if the decontamination is checked and judged as insufficient by us. It is your responsibility to bear the cost of sterilization or decontamination. Note that we cannot repair or inspect the rotor or the accessory unless sterilization or decontamination is completed.

9. Rotor Retirement

After many years of use, there will be inevitably some corrosion or stress corrosion. At some points, the combination of such damage and metal fatigue could make the rotor vulnerable to a failure.

Although a rotor may appear to be in a good condition, you should follow the rotor retirement recommendation shown below.

Rotor	Material	Retire After Years
Angle rotor	Titanium alloy	12
Vertical rotor	Aluminum alloy	10
Swing rotor	Titanium alloy Aluminum alloy	10
Zonal rotor	Titanium alloy	10
Continuous flow rotor	Titanium alloy	10

MEMO

MEMO

MEMO

Eppendorf Himac Technologies Co., Ltd.

1060, Takeda, Hitachinaka City
Ibaraki Pref., 312-8502 Japan

URL: <https://www.himac-science.com>

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Copyright © 2020 Eppendorf Himac Technologies Co., Ltd. All rights reserved.